MICROCLAVE

4001404

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GENERAL INFORMATION

- 1) *Handle the parcel with care. Unpack and check that the contents coincide with the packing-list. If any part is damaged or missing, please advise the distributor immediately.*
- 2) *Do not install or use the equipment without reading this handbook before.*
- 3) *This handbook must always be attached to the equipment and it must be available for all users.*
- 4) *If you have any doubts or enquiries, please contact with your supplier or J.P. Selecta's technical service.*
- 5) **IMPORTANT! J.P. SELECTA WILL NOT ACCEPT ANY APPARATUS TO BE REPAIRED IF IT IS NOT DULY CLEANED.**
- 6) *If any modification, elimination or lacking in maintenance of any device of the equipment by the user transgress the directive 89/655/CEE, the manufacturer is not responsible for the damage that can occur.*
- 7) *Do not use the apparatus with liquids which can give off vapours capable of making explosive mixtures.*

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CONTENTS

GENERAL INFORMATION
CONTENTS
PACKING LISTS
ACCESSORIES
TECHNICAL FEATURES
EQUIPMENT DESCRIPTION
INSTALLATION
OPERATION
MAINTENANCE
GUARANTEE
ELECTRICAL DIAGRAM
SPARE PARTS
"EC" CONFORMITY DECLARATION

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The standard equipment consist of the following components:

Descripción / Description	Código / Code	Cant. / Qty.
MICROCLAVE	4001404	1
Cestillo / Basket	1004770	1
Manual de instrucciones / Instruction Manual	80050	1

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* Baskets in perforated plate of AISI 316 stainless steel:

Código cestillo Basket code	Medidas cm. Dimensions cm	
	Ø	Altura Depth
1004770	21.5	15

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Voltage supply 115-230V 50/60 Hz according to the characteristics plate indications.

Código / Code		4001404
Capacidad litros / Capacity litres		7
Medidas útiles cm. Chamber dimensions cm.	Ø	24.5
	Alto / Depth	15
Medidas exteriores cm. Overall dimensions cm	Alto / Height	40
	Ancho / Length	28
	Fondo / Width	33
Consumo / Consumption W		1000
Peso / Weight Kg		13

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EQUIPMENT DESCRIPTION

For temperatures of 121°C and 131°C (1 and 1.7 Kg/cm²). Top and outer casing in AISI 304 stainless steel, chamber and lid in special aluminium alloy and a teflon-coated inner casing.

Built in a simple safety bayonet lock system and a safety valve.

INSTALLATION

Place the autoclave near a suitable electric mains.

Immobilise the autoclave with the stops provided in the front part of the apparatus.

Always use decalcified water.

Fix a hose in the water/steam exit nipple of the machine, and also fix the other end to a drain or deposit without **OBSTRUCTING THE FLOW**.

Do not install the autoclave in places where inflammable liquids are stored or in zones with special protection.

Before opening the cover **ALWAYS** check that there is **NOT PRESSURE**. The pressure gauge indicator installed in the control panel must read 0kg/cm².

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Be sure that the voltage supply is the same as the one indicated on the characteristics plate.

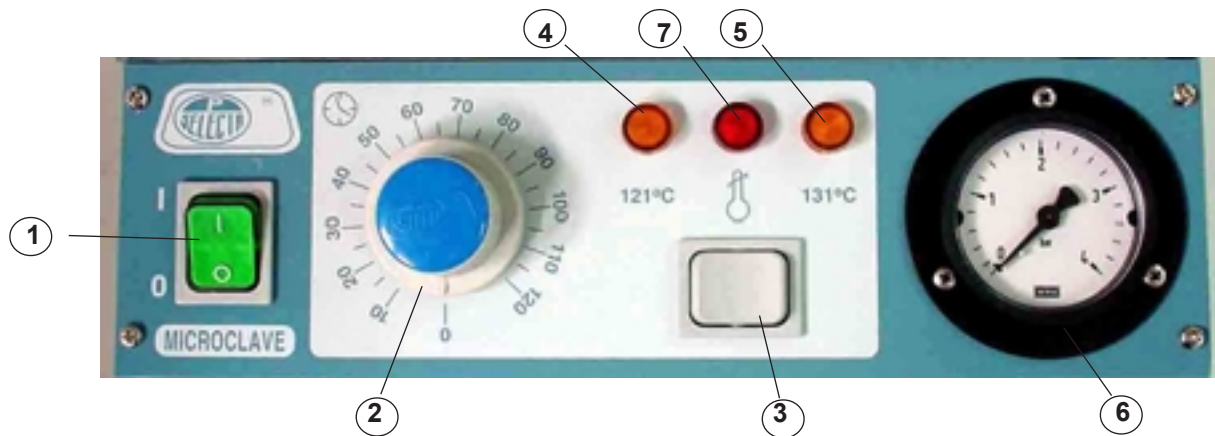
Do not use the autoclave if it is not earthed.

NEVER exceed the maximum pressure of 1.7 Kg/cm²

From time to time revise the functioning of the safety valve.

If you change the plug bear in mind the following:

Blue cable: Neutral.
Brown cable: Phase.
Yellow/green cable: Earth.

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OPERATION

STARTING UP:

1. Open the cover turning it anticlockwise. Fill the tank of the autoclave (better with decalcified water) up to the brim level (approx. 1 liter).
2. Place the material to be sterilized.
3. Place the side valve to the STEAM position.

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4. Put the cover, softly press down and turn clockwise until the two handles stay one above the other.

5. Seleccionar la temperatura de esterilización, a través del selector (3), teniendo presente la relación existente entre la temperatura y la presión de vapor saturado (Ver tabla 1).
5. *Select the sterilization temperature with the switch selector (3) bearing in mind the relation between temperature and pressure of the saturated steam. (See table 1).*

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11. To empty the inner water, place the handle to DRAINAGE position.

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SAFETY

* SAFETY THERMOSTAT:

When abnormal increase of the temperature or a considerable decrease in the water level occurs (with the consequent overheating of the heater element), the safety thermostat starts working blocking the heating of the machine and the ALARM indicator lamp (7) lights up. To deblock the machine proceed as follows:

1. Disconnect the autoclave from the mains.
2. Open the STEAM valve.
3. Let it cool.
4. Open and fill the tank with water up to the working level.
5. Reset the safety thermostat placed at the rear bottom part.
6. Start up the autoclave.

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If the problem continues, you must call the Technical Service.

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Said valve is gauged to initiate the opening at 2kg/cm², and only begins to work if the pressure switch fails.

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MAINTENANCE

Before removing the cover disconnect the apparatus from the mains.

The manipulation of the internal electronic circuits of the apparatus by unauthorized personnel can cause irreparable damage. Take it to one of the J.P.SELECTA authorized technical services.

CLEANING:

For the cleaning of the different parts use the following products:

Cleaning of stainless steel: Alcohol

Cleaning of plastic: Alcohol with cotton duster.

Cleaning of pipes with lime: Solution of Hydrochloric acid 50% and water 50%.

OTHER RECOMMENDATIONS:

Is advisable to change the water after 50 uses or once a month.

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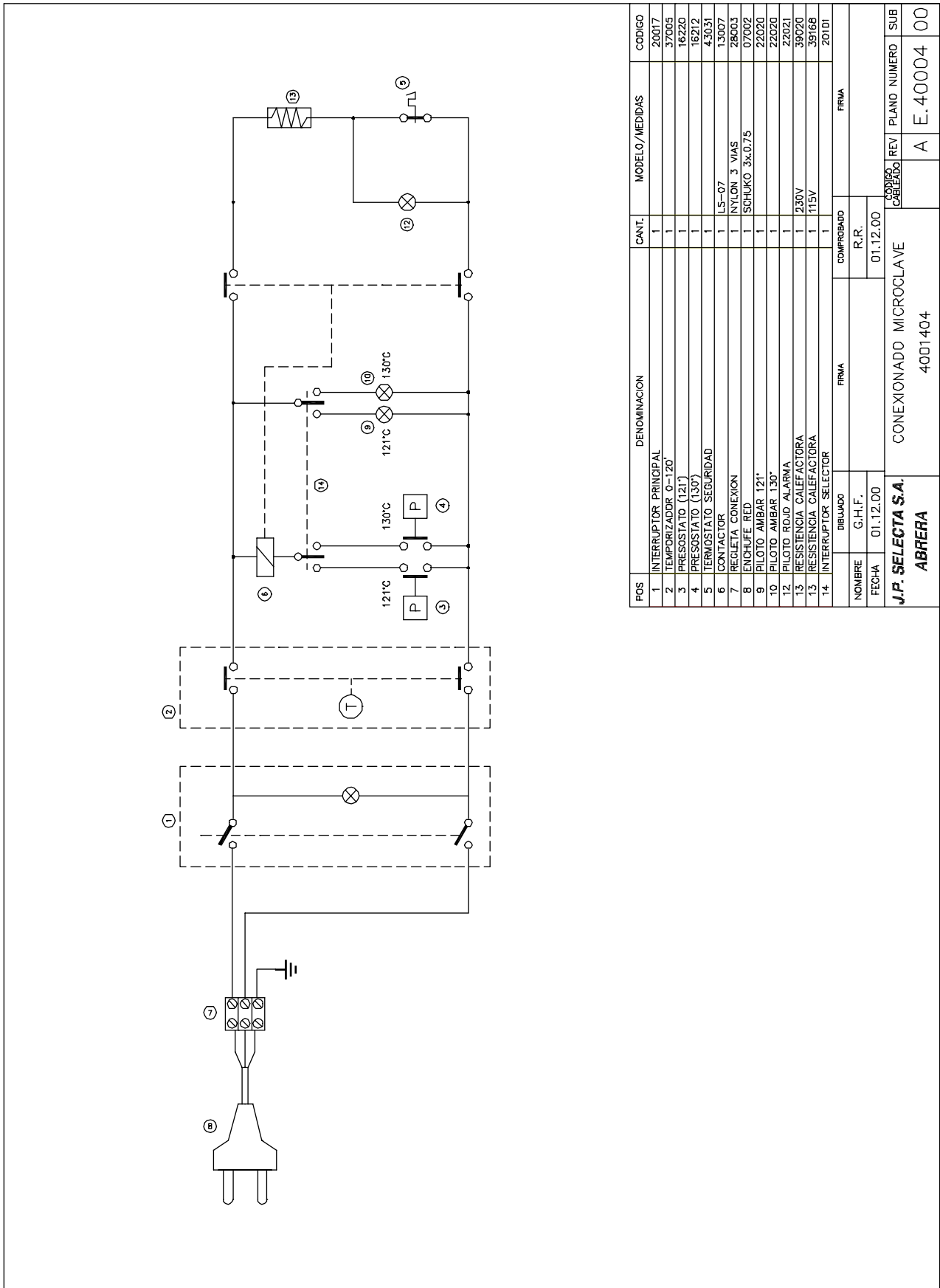
Annual revision of the whole function of the autoclave by qualified personnel is advisable.

GUARANTEE

This product is guaranteed for one year. The guarantee does not cover damage caused by incorrect use or causes beyond the control of J.P. SELECTA, S.A.

Any manipulation of the apparatus by personnel not authorized by J.P. SELECTA, S.A. automatically cancels the guarantee.

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To guarantee the safety of the equipment, the spare parts must be purchased from J.P.SELECTA, S.A.

Descripción / Description	4001404
Gradilla / Tube rack	5599
Mango tapa / Cover handle	15469
Manómetro / Manometer	16002
Presostato / Pressure switch	16220
Interruptor principal / Main switch	20017
Interruptor selector / Selector switch	20061
Junta tapa / Gasket cover	21017
Piloto ámbar / Ambar neon lamp	22020
Piloto rojo / Red neon lamp	22021
Válvula seguridad / Safety valve	34202
Reloj / Timer	37005
Resistencia / Heating element	39020

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